## KJM 5900 Oppgavesett 2:

# Gamma deteksjon og spektroskopi

[Exercise set 2: Gamma Detection and Spectroscopy]

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[This exercise is the first "mappeinnlevering" to KJM 5900. The submitted report count **10% on your final mark. The deadline is Friday 1<sup>st</sup> October at 4 pm. Failure to meet the dead line means you fail on this part of the exam.**]

The English translation of this text is located at the end, after the Norwegian text. If you do not want the translation you only need to print out the four first pages.

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## English translation

### How to perform and report the exercise

The report shall be written as a scientific paper. The paper should form a stand alone document, i.e. it should be possible to read it without having to refer to other sources to understand its content. However, it is quite alright, and also demanded, to make references to the course book or other sources when you make certain claims or draw conclusions which is not supported by your own work and results. As a guide to content and style, we have links to a few scientific papers on the main web-page.

The report should in most cases consist of the following parts:

- Title, author, date and publication place (UiO, Blinder).
- Abstract - write a short summary of your report.
- Introduction - Put your work in perspective. What is the aim of the exercise?
- Theory - Explain the background for the exercise (by this you answer part 2A and the theory part of 2B).
- Experiments and results - Describe what you have done and how (by this you answer the experimental part of 2B and 2C).
- Discussion and analysis - Describe how you analysed your measured data to reach your conclusions.
- Conclusion - What have you done and learned? Summarise your main results and draw conclusions.
- References - Here you list the sources you have used.

The report should not be longer than 5 typed A4 pages (with normal margins and 12 points font-size), exclusive figures and tables. Most likely your report will be better if it's shorter than this. A long length is not synonymous with quality....

Figures are nearly always of great help in explanations. But they can be quite time consuming to create. Therefore, you should carefully consider what kind of figures and how many you need. At the KJM 5900 homepage there is a pointer to a directory with a number of general figures. If you want to feel free to use this in your report. The directory is password protected: Username is *kjm5900* and the password is *curie*.

Read through this exercise description completely, including appendix A and B, before you start. It's also smart (necessary) to work through the guides on how to use the simulation program. You'll find links to the guides at the network page, as usual.

Submit the report as an attachment to e-mail. We prefer PDF formatted files, but if you use MS Word, LaTeX or WordPerfect then you can submit the file in these formats also. In that case we will generate a PDF file for you and send it back to you as a check.

### Introduction to this exercise

Absorption of  $\gamma$  radiation occur by three different effects: The photoelectric effect, the Compton effect and pair formation. These effects are described in the course book in chapter 6.4.

A NaI crystal will emit light when  $\gamma$  radiation is absorbed. If we optically couple the crystal to a sensitive light-detector (a photo multiplier tube), we can measure the amount of light emitted and therefore also the  $\gamma$  radiation. In this way we can construct a  $\gamma$  detector and we call it a NaI-scintillator detector (to scintillate means emits light). Such detectors are described in the course book in chapter 7.3.

The amount of light is proportional to the amount of energy deposited in the detector. The light detector is so fast that it's able to measure the light from each individually absorbed  $\gamma$  ray. By analysing

the pulse height from the detector (which is proportional to the amount of light) we can make a picture of the distribution of energy deposited in the detector. I.e. we make a histogram where each column represent the number of events within a small energy range - we will get a plot of event intensity as a function of energy deposited in the detector. Such a histogram we call a *spectrum*.

If we have a source which only emits  $\gamma$  rays with a single energy, its spectrum should look like Figure 1. However, if we measure such a source with a NaI scintillation detector, the spectrum we get look like Figure 2. The reason is of course the manner  $\gamma$  rays are absorbed in the detector. A large fraction of the  $\gamma$ s will be absorbed by the Compton effect and this severely distorts the spectrum.

In addition, the detector do not measure the amount of energy deposited with 100% accuracy. The error in the measurement will widen the peaks and generally smooth the spectrum. Thus, details will be more difficult to distinguish. The detector error is generally discussed in terms of the detectors *response function*.

## Exercise 2A: $\gamma$ emission from radio nuclei

Describe what  $\gamma$  radiation is and how it interacts with matter. Describe how a NaI crystal and a photo multiplier tube (light detector) can be used for measuring  $\gamma$  radiation.

## Exercise 2B: Detection of $\gamma$ radiation

In a real detector one can not turn the different absorption effects on and off - they will always be on. However, in this exercise we will use a program which simulated a real detector and then you have full control of which effects should be active or not. You can even vary the intensity of the different effects. The detector-simulation application you can download from the KJM 5900 web-page and run it as a stand-alone application. The link to the application is called "[Gamma Detector Simulation](#)". *Make certain that the computer you use has the LabView run-time library installed, otherwise the application will not run.* This and other practical things are explained in the instructions, which you access from the web-pages.

The application enables you to control four detector properties: The response function (responsible for the measurement error) and the absorption coefficients for the photo electrical effect, the Compton effect and pair formation. Furthermore the application let you chose if your source should emit  $\gamma$  rays with one, two or three different energies. And you can vary the intensity of which each energy is emitted with.

Perform the following experiments with the program, where you select only a single  $\gamma$  energy with 662 keV of energy:

- Make a "perfect" spectrum, as it would be measured if only the photo electrical effect was present and no detector error. I.e. set the application's response function to 0, the photo electrical effect to 100%, the Compton effect to 0% and pair formation to 0%.
- Make a spectrum which has no detector error (response function = 0), but only absorbs  $\gamma$ -rays with the Compton effect (photo electrical effect = 0%, Compton effect = 100% and pair formation = 0%). Compare this spectrum to the one you made in a). Which one is most suitable for analysing the radiation energy and intensity from a  $\gamma$  source?
- Make a spectrum where only the photo electrical effect is at play, but select a response function of 100. Can you explain the shape of the response function? Tip: Some times the primary electron (the one which the  $\gamma$  ray transferred its energy to) will not excite the NaI crystal as much as it should have done. This

might be due to impurities in the crystal or that the electron exits the crystal before it has been stopped completely. In other words we will have some events which can have anything from 0% and up to 100% of the correct energy. For example, an electron which deposit 50% of its energy and then escapes the crystal will be registered as a count with only half of the "correct" energy.

d) Make a spectrum with the response function set to 100, the photo electrical effect equal to 50% and the Compton effect equal to 50%. Compare this to a spectrum from a real NaI detector (download it from the KJM 5900 web page).

For each experiment you should write down the parameters you have used and store the spectrum as an image file. In this way you can incorporate your spectra in your report. You store the spectra by right clicking it and then choosing the menu command "Export Simplified Image". It's the same procedure as when you pick pictures from web-pages in most browsers.

### **Exercise 2C: Analysing spectra with multiple $\gamma$ -lines**

Use the simulation application you used in 2B, but this time you select three different  $\gamma$  ray energies: 400 keV (10% intensity), 450 keV (50%) and 800 keV (100%).

a) Make a spectrum with an ideal detector (response function = 0, photo electrical effect = 100%, Compton effect = 0% and pair formation = 0%).

b) Compare this spectrum with what you get if the response function is 10 and 100.

c) What happens if you turn on the Compton effect?

### **Conclusion**

In your report's conclusion, you should try to explain what you have learned about  $\gamma$  spectroscopy. E.g. how does the detectors energy resolution (the response function) influence the quality/uncertainty of the spectrum? And what would be the advantage of selecting a detector made in such a way that the photo electrical effect is as dominant as possible?

## Appendix A: Help!

Nobody expect you to be an expert on  $\gamma$  spectroscopy. Try to write a report which demonstrates that you have understood the main issues concerning detection of  $\gamma$  rays. In particular it should convince the reader that you understand how  $\gamma$  rays are absorbed, how a NaI detector works and how you should interpret a  $\gamma$  spectrum. If you at the same time are able to express yourself in a precise manner, avoid too much "spoken language" in your text, have a good disposition and that your report "stand by itself", i.e. that it can be read as a stand-alone text, then you most likely will deliver a good report.

Do not refer to this exercise text as the reason for doing certain things - you should explain the reason for doing it, not claim that your teachers forced you to do it... Understanding *why* you are asked to perform the different operations is a large part of learning about the subject!

It might be smart to obtain further information from other sources than the course book. For example, you could make some searches on the net with key words like "Compton edge gamma spectroscopy" or something similar. Most likely you will find something useful. Beware of getting into too much detail, though - first of all we want you grasp a basic understanding of  $\gamma$  spectroscopy. Keep in mind that there is a page limit!

In the counting room in the course lab we have a small collection of text books. Feel free to pop in and browse through these books (but do not take them with you!).

Keep an eye on how much time you use for this exercise. Certainly you have other courses and tasks which demand attention and even if the report is 10% of your final grade, it's not so important that you should ignore everything else!

If you have problems with the simulation application you should notify the supervisors immediately. If you are at Blindern the counting room at the course lab has PCs which has the LabView run-time installed. You can use these PCs for this exercise (or any other KJM 5900 related tasks) as much as you want.

## Appendix B: Criteria used for the evaluation

1. Do the report meet the demands of scientific writing?
  - a. Disposition and selection of content.
  - b. Language (should not be too oral).
  - c. Use of sources.
2. Readability and completeness
  - a. Do the student put forward her/his arguments in a manner which is easily understood?
  - b. Is the amount of theory well balanced to the experimental part? (The theory should only support the experimental part, not be a teaching book.)
  - c. The correct use of figures and tables - they should aid in the description and organisations of the data and analysis presented. They should form a natural part of the report.
3. Has the student answered the questions asked?
  - a. Exercise 2A.
  - b. Exercise 2B.
  - c. Exercise 2C.
  - d. Is the students conclusions correct?

Each criteria listed above counts with equal weight in the final evaluation.